



Directions for Digital Microscopes
USB Shot
SAVING

1. Create a folder on your desktop by clicking on the desktop, then go to File, and New Folder

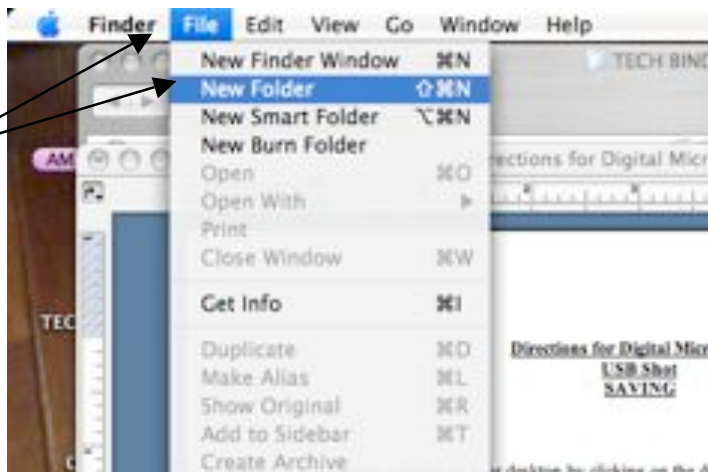


2. When the new folder appears on your desktop, type and label it microscope.
(if you type and nothing changes. Click on the label until you see a cursor, then type)

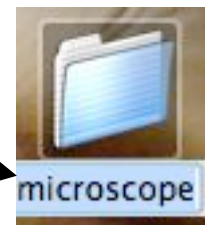


3. With the microscope plugged in, open USB SHOT.

4. Make sure that you are in Snap Shot Mode. If you are not sure click on Snap Shot.

5. Click on the folder icon.

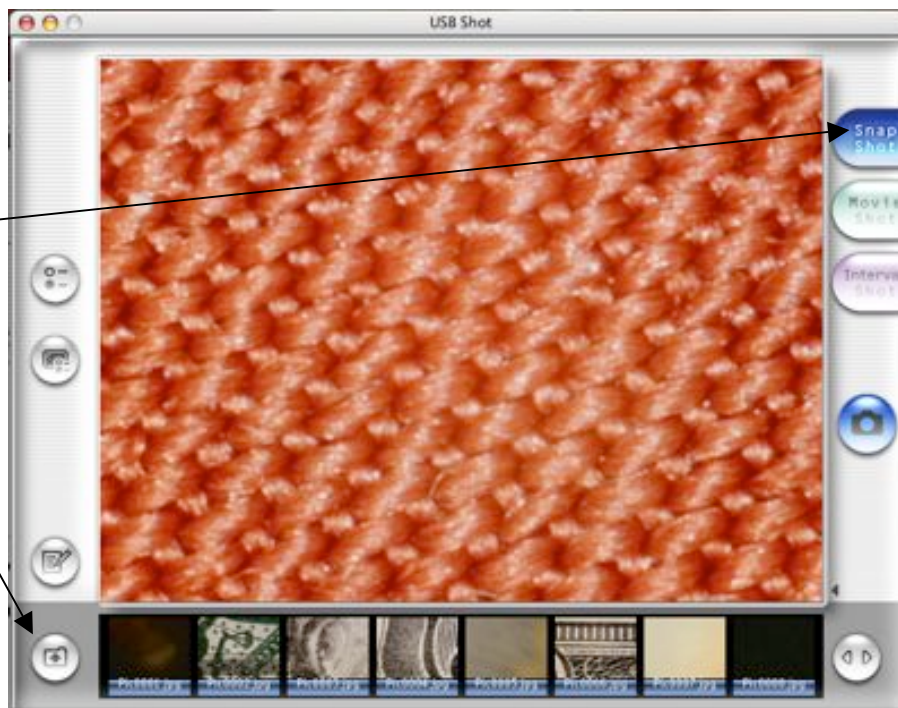

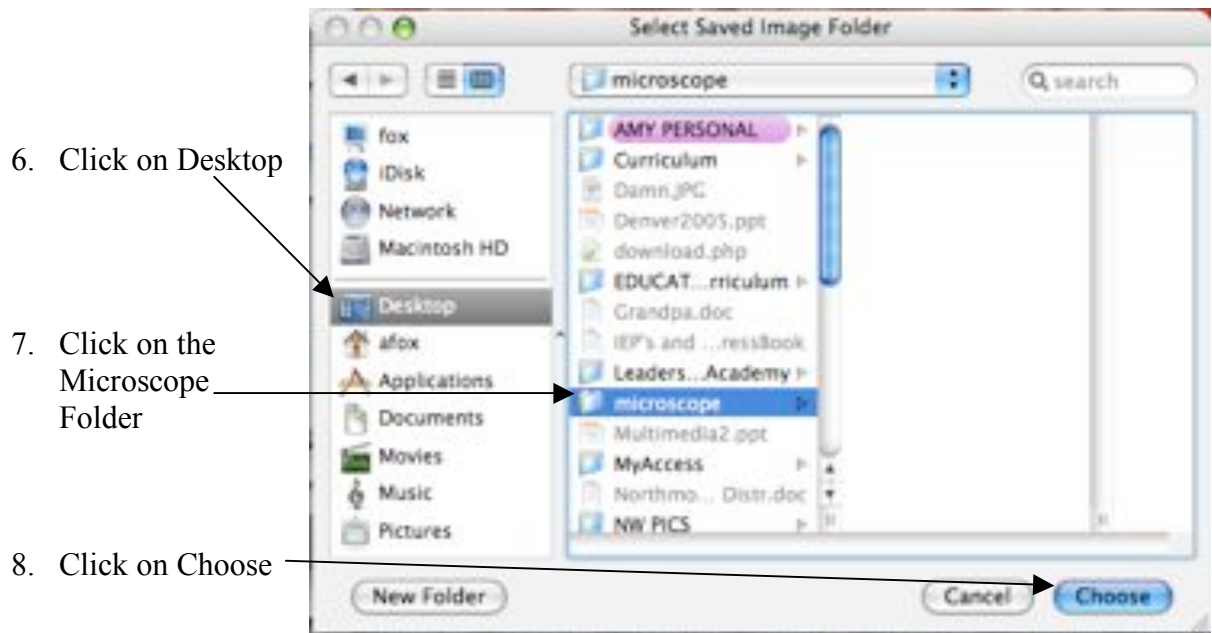

Now your snap shots will be saved to this folder.

***** Repeat step 3 to 8 for Movie Shot and Interval Shot as well.*****